

Multi-class pesticides analysis in challenging vegetable matrices using fast 5 msec MRM with 15 msec polarity switching

ASMS 2012 WP27-575

Yuka Fujito¹, Yusuke Inohana², Kiyomi Arakawa², Ichiro Hirano² ¹Shimadzu Analytical & Measuring Center, Inc., Kyoto, Japan ; ²Shimadzu Corporation, Kyoto, Japan Multi-class pesticides analysis in challenging vegetable matrices using fast 5 msec MRM with 15 msec polarity switching

Introduction


Many regulatory authorities have established multi-class residual pesticides methods for the analysis of vegetables, fruits and other food stuffs. There is, however no global agreement on the provision of a target list of pesticides and this presents a risk with products moving between different regulatory requirements. In order to eliminate this risk, food safety laboratories need to ideally screen as many compounds as possible in a single run which may reach maximum residual limits (MRL); typically 10 ppb in food matrices.

In this study, we report the application of ultra-fast 5 msec MRM with 15 msec polarity switching for the analysis of 138 pesticides in vegetable matrices (72 and 66 compounds measured by LC-QqQ and GC-QqQ in the European Union Reference Laboratory (EURL) method).

Materials and Methods Sample Preparation (QuEChERS EU method)



Food sample

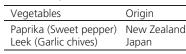

Step 1 : Sample Extraction

1. Homoginize vegetables with food processor and homoginizer

2. Weigh 10 g homoginized sample



4. Centrifuge for 5 min. at 4000 rpm (Extract 1)

*1 : Citrate Extraction Tube (SIGMA ALDRICH)

Step 2 : Sample Clean up



*2 : PSA/ENVI-Carb SPE Clean Up Tube 2 (SIGMA ALDRICH)

LC/MS/MS analysis

Analytical Conditions

Column	: Shim-pack XR-ODSII (75 mm x 2 mml.D., 2.2 um)
Mobile phase	: A ; 2 mM ammonium formate containing 0.1%
	formic acid – water
	B ; Methanol
Gradient program	: 5% B (0-2.5 min.)→55% B (2.51-6 min.)→80%
	B (6.01-12 min.) →100% (12-15 min.)→5%
	(15.01-20 min.)
Flow rate	: 0.2 mL / min.
Column temperatur	re : 40°C

MS : LCMS-8040 Triple quadrupole mass spectrometer

Ionization	: ESI (Positive / Negative)
lon spray voltage	: +4.5 kV / -3.5 kV
MRM	: 276 MRM transitions (2 MRMs / compound)
	Dwell time 5 msec. / Pause time 1 msec.

Features of LCMS-8040

- 5 times higher sensitivity compared to LCMS-8030
- An ultra fast scan speed of 15000 u / sec.
- An ultra fast polarity switching of 15 msec.
- An ultra fast MRM transition speed of 555 ch./ sec.



Fig. 1 LCMS-8040 Triple Quadrupole Mass Spectrometer

Multi-class pesticides analysis in challenging vegetable matrices using fast 5 msec MRM with 15 msec polarity switching

Results MRM of pesticide standards

Table 1	100s of 138	pesticides in the FUR	L method by LCMS-8040
Tuble I	LOQJ 01 130	pesticides in the Lon	

Technique (on the EURL method)	LOQs < 10 ppb	LOQs > 10 ppb	Not Ionization
by LC-QqQ	72 (100%)	0 (0%)	0 (0%)
by GC-QqQ	47 (71%)	6 (9%)	13 (20%)

• 71 % of pesticides measured by GC-QqQ in the EURL method could achieved excellent LOQs (0.08-10 ppb) .


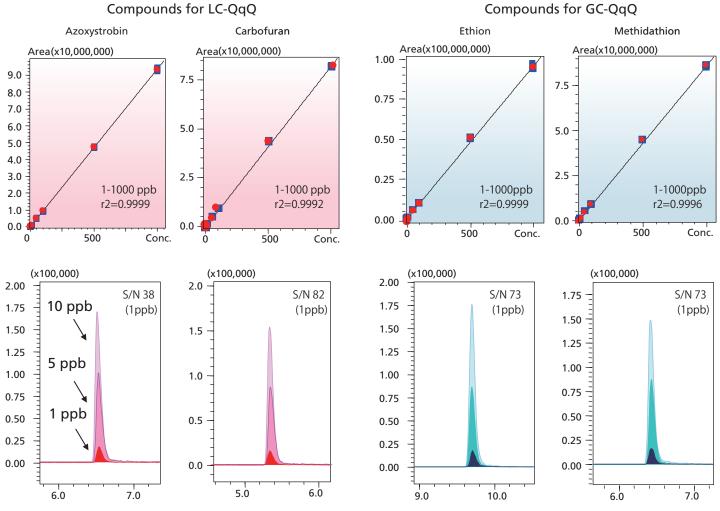


Fig. 2 Calibration curves and MRM chromatograms of typical pesticides

Multi-class pesticides analysis in challenging vegetable matrices using fast 5 msec MRM with 15 msec polarity switching

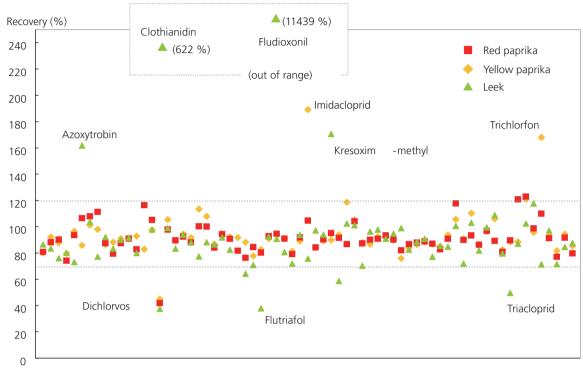

Recovery of pesticides in vegetable matrices

Fig. 3 Recovery of pesticides in the vegetable matrices (5 ppb spiked)

- Approximately 90% of pesticides represented good recoveries in the range of 70-120% in all studied matrices.
 Some pesticides indicating up to 120% of recovery were also detected in the matrix blank (Fig. 4).
 - ----- 5 ppb standards spiked ------ Matrix blank

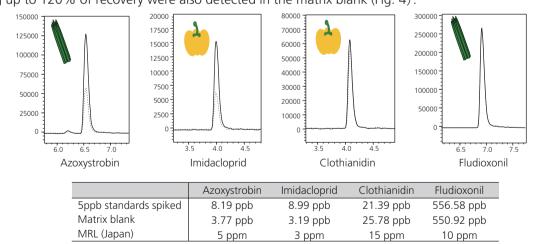


Fig. 4 MRM Chromatograms and quantitative results of 4 pesticides

Conclusion

- MRL of pesticide in vegetable matrices pre-treatment by QuEChERS method could be measured successfully using fast 5 msec MRM with 15 msec polarity switching.
- Majority of pesticides being suggested to GC-QqQ technology on the EURL method was successfully covered by LC-QqQ.



Shimadzu Corporation www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures. The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

© Shimadzu Corporation, 2012